CAPSULE FILLING OPERATION - STEPS



Step 1

Load the capsule body, Put the body plate in a flat.



Step 2

Put the frame of the arrangement plate on the body plate.



Step 3

Put about 100 Capsules in the frame.



Step 4 pick up the plate body and arranged wheel swinging around.



Step 5 Capsule body will fall into the plate hole one by one, then, pour the extra capsule body out.



Step 6 Take the wheel arrangement plate, put the capsule body installed on the board aside.



Step 7

Filling power in the capsule body.



Step 8

Scrape the board back and forth on the body, When the capsule body is full on powder.



Step 9

Use powder pressing plate, filling capsules.



Step 10

Scraping plate excess powder.



Step 11

Turned overthe the cap plate and middle plate to cover the body plate.



Step 12

Lock the capsule body and the capsule cap.



Step 13

Remove the cap plate.



Step 14

Remove the middle plate, pour out the capsule from the locking middle plate.